

# AAA-3500


## Automatic Analyzer of Amino Acids



- ▶▶ AAA 3500 is a compact device designed for determination of amino acids and biogenic amines. Thus, it can be used in the fields of biochemical research, research of human and animal nutrition, medical diagnostics, control of drugs, food and feedstock. The robust and sensitive method of ion exchange chromatography with post-column ninhydrin derivatization provides high accuracy and reproducibility of analyses. AAA 3500 will provide results with maximum accuracy and reproducibility in continuous operation, saving purchasing, operation and maintenance costs.

## Compact device

AAA 3500 is a compact device with a unified control panel, which increases user convenience and facilitates control and handling of the device. AAA 3500 is normally controlled by a computer, the user may control the settings of the device with the use of a touch display as well.

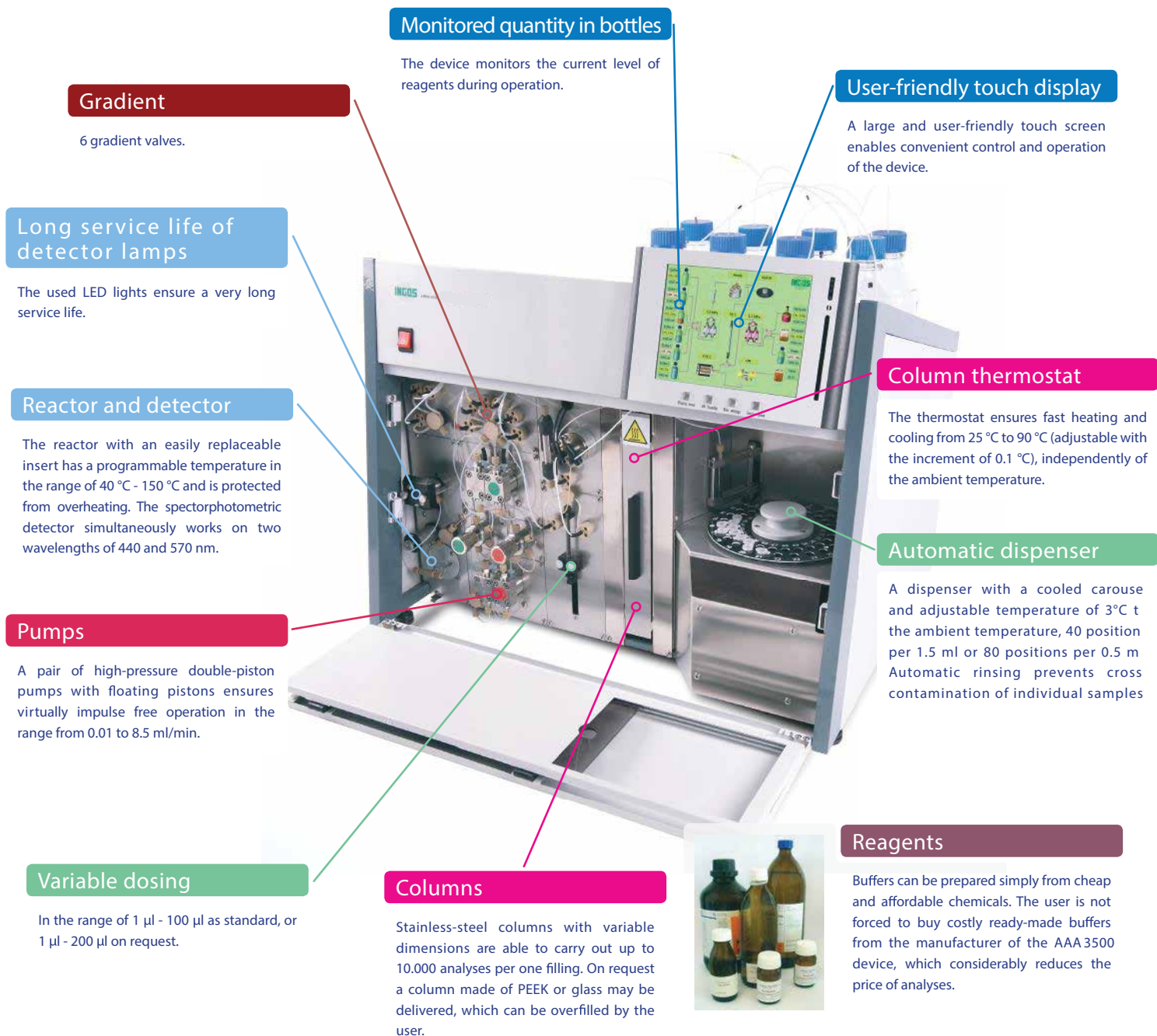


## Robust and sensitive method

AAA 3500 uses the principle of ion exchange chromatography with post-column NHD derivatization. Using this method you can safely determine amino acids in hydrolysates of proteins, peptides as well as free amino acids in physiological solutions and extracts, biogenic amines and some antibiotics.

## Saves on costs

The user can minimize the operation costs by producing reagents with the use of simple software from cheap and affordable chemicals (there is no need to buy overpriced ready-made buffers).

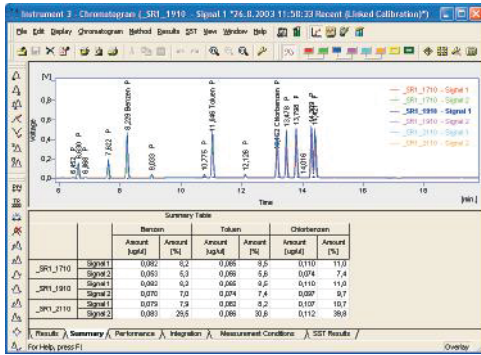

## Clarity control software


AAA 3500 is controlled by the Clarity software of the DataApex Company. This software ensures convenient and intuitive operation of the devices as well as work with data and their export to various formats. Clarity makes it possible to connect other equipment to AAA 3500, e.g. a fluorescence detector.

The device can be controlled remotely via the Internet or Maintenance diagnostic activities can also be carried out remotely.


## High accuracy and reproducibility of analyses

AAA 3500 provides analyses with a low level of noise and drift. The ATL Company is able to produce a shortened or modified program for virtually all ninhydrin positive substances on the customer's request.


### Determination of free amino acids (standard)

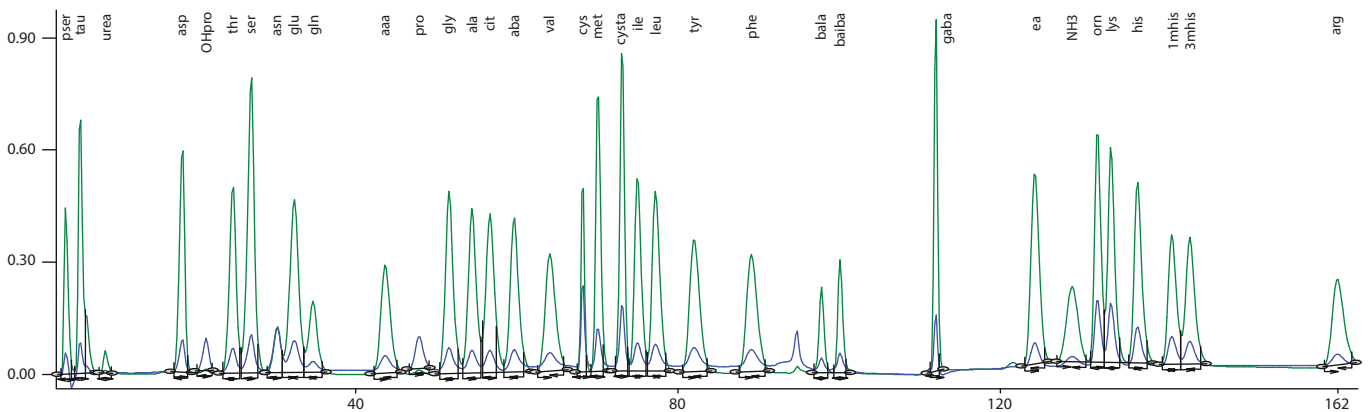


### Hydrolysate analysis (standard)

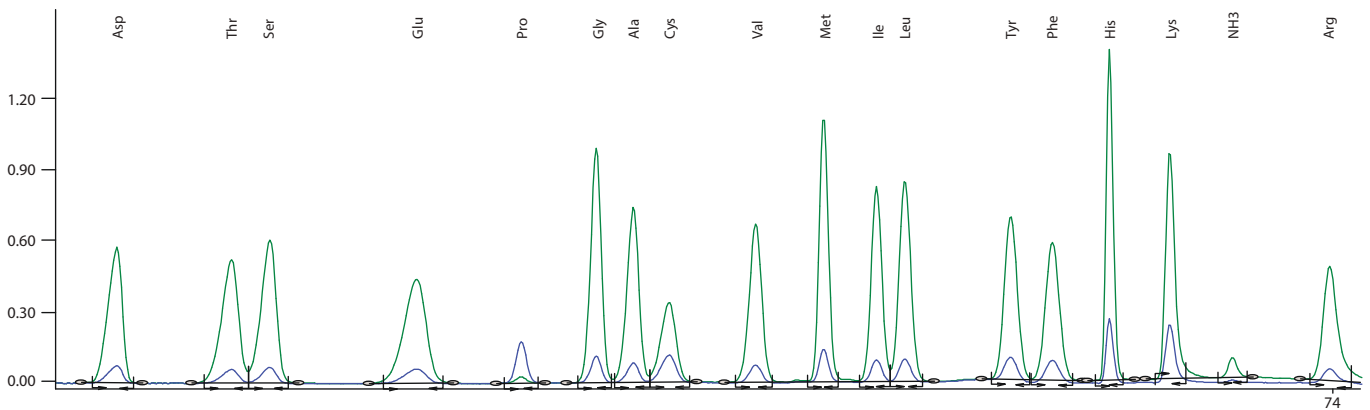

### Shortened determination of sulphurous amino acids (cystine and methionine)

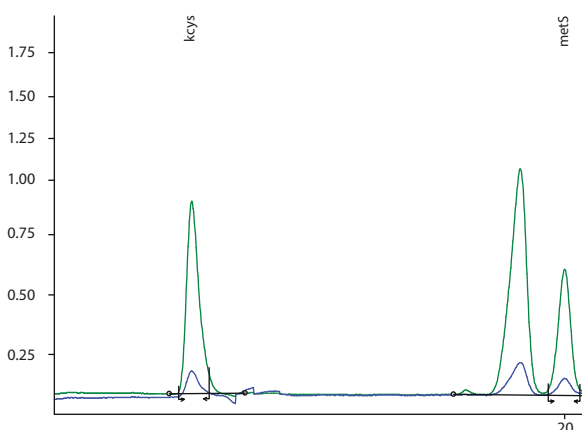

### Determination of biogenic amines

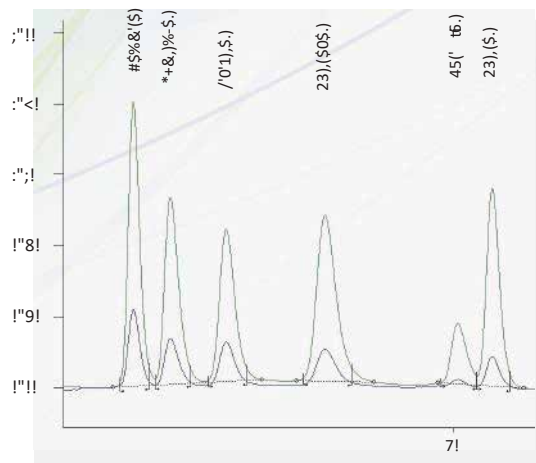

## Technical specifications

### Automatic Analyzer of Amino Acids 3500

Noise	1.7 x 10 AU/s
Drift	2.0 x 10 <sup>-5</sup> AU/hod
Reproducibility at 10 nmol	Retention times: 0.3 % (arg) Peak area: 1 % (ser, gly, his)
Pumping system	High pressure impulse-free pumps (pair of floating pistons)
- flow	0.01 to ≤10 ml/min
- pressure	0 – 25 MPa
Dispenser with cooling	
- no. of samples, cassette A	40 x 1.5 ml
- no. of samples, cassette B	80 x 0.5 ml
- variable dosing	1 - 200 µl with 1 µl increment
Cooling	Peltier
Temperature	Adjustable 3 °C (ambient temperature)
Columns	
- dimensions	4 x 250 mm or 4 x 150 mm
- material	stainless steel, PEEK, glass
Column thermostat	
- temperature	25 °C – 90 °C
- cooling and heating	Peltier
- setting accuracy	0.1 °C

Degasser	6-channel degasser
Double-channel photometer	400 and 600 nm
Cuvette volume	5 µl
Reactor	40 – 150 °C (overheating protection)
PC communication	LAN
Processing of results	Clarity
Colour display	8" diagonal
Power input	230V, 50Hz
- start	280 VA
- operation	120 VA
- standby	70 VA
Dimensions (W x H x D)	710 x 600 x 560 mm
Weight without bottles	46 Kg
Operation time without the user's intervention	approx. 7 days
Other applications	Biogenic amines Sugar analysis for reduced sugars

The ATL Company offers  
e.g. these laboratory instruments:





## HPLC Servicing, Validation, Trainings and Preventive Maintenance :

**HPLC Servicing** :HPLC Servicing : We have team of service engineers who can attend to any make of HPLC promptly @the most affordable cost.

**Trainings** :We also take up preventive Maintenance to reduce downtime of HPLC's Trainings.

**AMC's/CMC** :AMC's/CMC :We offer user training both in-House and at customer sites on HPLC principles, operations, trouble-shooting.

**Validations** :Validations :We have protocols for carrying out periodic Validations as per GLP/GMP/USFDA norms.

**Instruments** :Instruments :We offer instruments/Renting Services Modules like pumps,detector etc. on Rent.



## About Analytical Technologies

Analytical Technologies is synonymous for offering technologies for doing analysis and is the Fastest Growing Global Brand having presence in at least 96 countries across the global. Analytical Technologies Limited is an ISO:9001 Certified Company engaged in Designing, Manufacturing, Marketing & providing Services for the Analytical, Chromatography, Spectroscopy, Bio Technology, Bio Medical, Clinical Diagnostics, Material Science & General Laboratory Instrumentation. Analytical Technologies, India has across the Country operations with at least 4 Regional Offices, 6 Branch Offices & Service Centers. Distributors & Channel partners worldwide.

## Our Products & Technologies



UV/VIS  
Spectro 2080+  
Double Beam



Infra FTIR



Optima Gas  
Chromatograph  
3007



Optima Gas  
Chromatograph  
2979 Plus



Flash  
Chromatograph



Atomic Absorption  
Spectrophotometer



Liquid Particle  
Counter



Optical Emission  
Spectrophotometer



DSC/TGA



Semi Auto Bio  
Chemistry Analyzer



HEMA 2062  
Hematology  
Analyzer



Micro Plate  
Reader/Washer



URINOVA 2800  
Urine Analyzer



Total Organic  
Carbon 3800



Fully Automated  
CLIA



NOVA-2100  
Chemistry Analyzer



PCR/Gradient PCR/  
RTPCR



TOC  
Analyzer



Laser Particle  
Size Analyzer



Ion Chromatograph



Water purification  
system

## Regulatory compliances



## Corporate Social Responsibility

Analytical Foundation is a nonprofit organization (NGO) found for the purpose of:



# Analytical Foundation

1. Research & Innovation Scientist's awards/QC Professional Award : Quality life is possible by innovation only and the innovation is possible by research only, hence ANALYTICAL FOUNDATION is committed to identify such personalities for their contributions across various field of Science and Technology and awarding them yearly. To participate for award, send us your details of research / testing / publication at [Info@analyticalfoundation.org](mailto:Info@analyticalfoundation.org)

2. Improving quality of life by offering YOGA Training courses, Work shops/Seminars etc.

3. ANALYTICAL FOUNDATION aims to DETOXYFY human minds,souls and body by means of yoga, Meditation, Ayurveda, Health Care, Awards, Media, Events, Camps etc.

## Reach us @



HPLC Solutions   MultipleLabs   Analytical Bio-Med   Analytical Distributors   Analytical Foundation (Trust)

**Corporate & Regd. Office:**  
Analytical House, # E67 & E68,  
Ravi Park, Vasna Road, Baroda,  
Gujarat 390 015. INDIA

T: +91 265 2253620  
+91 265 2252839  
+91 265 2252370  
F: +91 265 2254395

E: [info@hplctechnologies.com](mailto:info@hplctechnologies.com)  
[info@multiplelabs.com](mailto:info@multiplelabs.com)  
[info@analyticalgroup.net](mailto:info@analyticalgroup.net)  
[info@analyticalbiomed.com](mailto:info@analyticalbiomed.com)

W. [www.analycalgroup.net](http://www.analycalgroup.net)  
[www.hplctechnologies.com](http://www.hplctechnologies.com)  
[www.multiplelabs.com](http://www.multiplelabs.com)  
[www.ais-india.com](http://www.ais-india.com)

Sales & Support Offices:  
across the country :  
Distributors & Channel  
partners World Wide

Note : Company reserves rights to add/delete/modify the contents / technical specifications of the catalogue without prior notice.