

AAS- 3810i

Atomic Absorption Spectrometer



EPCC / PRODUCTS / APPLICATION / SOFTWARE / ACCESSORIES / CONSUMABLES / SERVICES

Analytical Technologies Limited

An ISO 9001 Certified Company

www.analyticalgroup.net

AAS-3810i is the fifth-generation atomic absorption spectrometer developed. It combines the quality and precision that is known for with important new features. The AAS-3810i graphite furnace features a unique design with transverse heating, longitudinal Zeeman background correction and variable magnetic field intensity. The Zeeman background correction is suitable for high precision analysis of samples with complex interference matrices. Variable magnetic field intensity allows the operator to optimize field strength for each element, maximizing the Zeeman effect for accurate results every time.

The new standard for high performance Zeeman GF analysis available in a flame tandem configuration:



AAS-3810i atomic absorption spectrometer is a fully automated instrument with three main configurations: AAS flame analyzer, AAS Zeeman graphite furnace analyzer, and AAS-3810i flame and Zeeman graphite furnace tandem analyzer. Key features include:

- Transverse graphite furnace heating technology improves sample atomization efficiency and uniformity.
- Two background correction methods are available: longitudinal Zeeman effect correction and deuterium lamp correction. The operator can select the most suitable mode based on the sample.
- Unique adjustable magnetic field strength (0.6 - 1.1 Tesla, in 0.1 Tesla increments) allows for element specific optimization to achieve high sensitivity and precision.
- 8-lamp automated lamp turret with auto-alignment for quick element switching and lamp position optimization
- Code lamp recognition technology, compatible with both hollow cathode lamps and super lamps
- High efficiency super lamp power supply, newly designed for higher sensitivity while consuming less power and generating less heat

Unique functional design, easy to operate : Automated burner head height control

The burner head height can be automatically adjusted using a software-controlled motor. This feature is used to ensure that the burner head can be accurately and reproducibly positioned on the optical path to optimize the performance of the instrument.

The position of the burner head is stored with the analysis method parameters of each element in a multi-element flame analysis. During a sample run, each element will be analyzed at the optimal height without operator intervention. The following figure shows the sensitivity obtained when the burner head height of each element is separately optimized (purple) compared to the sensitivity obtained when the burner head height is optimized for only a single element, copper (cyan).

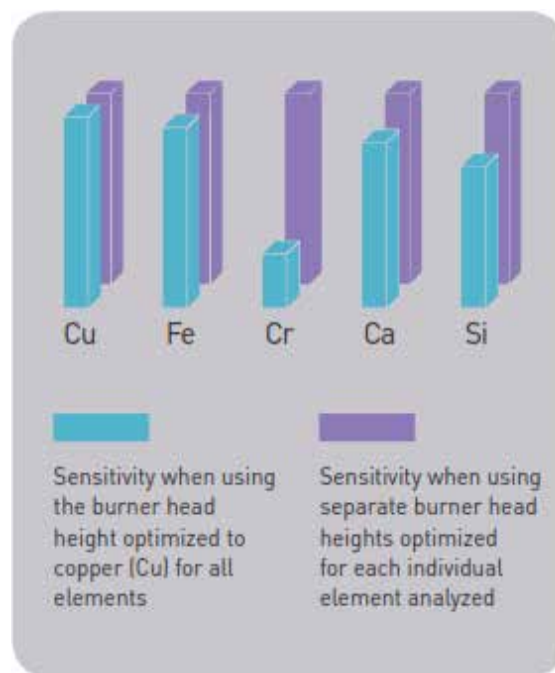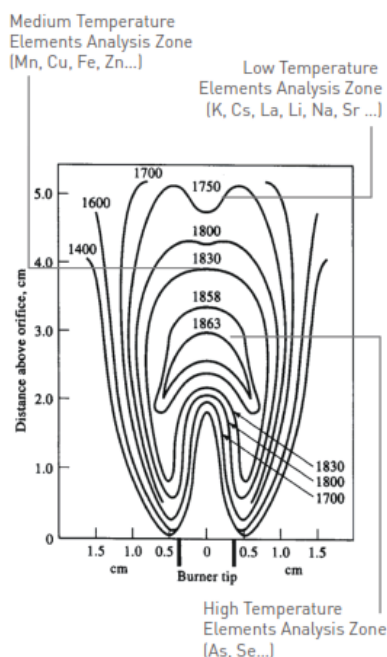

Super lamp power supply:

Super lamps are specially designed ultra-high intensity hollow cathode lamp that provide more optical energy. For certain elements, such as As, Se, Cd, Ni and Pb, super lamps can significantly improve the limit of detection, sensitivity and linearity. AAS 3810i atomic absorption spectrometer can be equipped with up to four super-lamp power supplies.

Coded lamp recognition:

By simply inserting the element lamp into the lamp holder, the software will automatically identify the element and position of the lamp. This feature completely eliminates the possibility of operator error when inputting lamp information into the element lamp list.

Graphite Furnace Viewing System:

The Graphite Furnace Viewing System uses a camera to view the flame or graphite furnace. This feature is indispensable for the development of graphite furnace methods. The operator can observe the entire analysis process from injection to atomization in real time. Using visual information, the desolvation, drying and ashing parameters can be set correctly to produce reproducible and accurate results.

Graphite furnace gas saving mode:

The intelligent control of the protection gas switch maximizes the effective use of protection gas and reduces waste when the gas is not needed, cutting down on the cost of operation.

Excellent performance, accurate results guaranteed

Cutting edge transverse heating, longitudinal Zeeman effect graphite furnace technology:

Compared with transverse Zeeman effect graphite furnace designs, the longitudinal Zeeman effect graphite furnace has the advantage of not needing a polarizer located along the optical path. This results in a two-fold increase in the amount of energy reaching the detector, which provides higher sensitivity and better correction linearity. Transverse graphite furnace heating has the advantage of higher temperature and greater temperature distribution uniformity, which promotes atomization of the sample with high efficiency and uniformity. Efficient graphite furnace design and precise temperature control provides the user with a reliable and precisely controlled analysis environment. This allows the AAS 3810i to perform automated multi-element, maximum heating rate guarantees optimal atomization conditions for all elements.

Unique Variable Magnetic Field Strength:

A constant magnetic field strength for all elements leads to severe limitations in analytical performance for many sample types. The AAS 3810i achieves optimal background correction and sensitivity by allowing optimization of the magnetic field strength for each element/matrix combination. The applied magnetic field can be varied from a low of 0.6 to a high of 1.1 Tesla in increments of 0.1 Tesla. This allows the operator to set the optimum field strength for each element, maximizing the Zeeman Effect as measured through the magnetic sensitivity ratio (MSR). Variable background correction allows the field strength to be adjusted to maximize both precision and sensitivity. This produces accurate results every time.

Unique double gas route:

The AAS 3810i features a unique double gas route design, with an inner gas route and an auxiliary gas route. Auxiliary gas is controlled through software. The addition of auxiliary gas promotes sample ashing and protects the graphite tube, allowing for more firings before replacement.

Two background correction modes:

The operator can select between Zeeman effect, and deuterium lamp background correction modes based on the requirements of a particular sample matrix.

Accessories:

AAS 3810i works with furnace autosampler, and flame/furnace autosampler. Also it can be used with hydride generator and other accessories.

Technical Specifications:

Optics:

Wavelength Range: 190 nm ~ 900 nm

Bandwidth: 0.1, 0.2, 0.4, 1.0 and 2.0 nm, automatic switch

Czerny-Turner Grating Monochromator

Wavelength Accuracy / Indication Error: ± 0.1 nm

Wavelength Repeatability: ≤ 0.05 nm

Grating: 1800 lines / mm

Blaze Wavelength: 250 nm

Baseline Stability: Static baseline stability ≤ 0.002 A / 30 min, dynamic baseline stability ≤ 0.003 A / 30 min

Resolution: Better than 0.1 nm

Graphite Furnace System:

Characteristic Concentration of Cd: $\leq 0.5 \times 10^{-12}$ g

Detection Limit: $\leq 1 \times 10^{-12}$ g

RSD $\leq 3.0\%$

Temperature Control: Max 20 step temperature program.

3 modes of temperature rise: step, slope and flat.

Heating Mode: power heating

Heating Rate: $\geq 2500^\circ\text{C} / \text{s}$

Flame System:

Characteristic Concentration of Cu: ≤ 0.02 $\mu\text{g} / \text{mL} / 1\%$

Detection Limit: ≤ 0.003 $\mu\text{g} / \text{ml}$

RSD $\leq 0.45\%$

Burner Head: Switchable 50 mm and 100 mm titanium burner head, titanium nebulizer and high performance glass nebulizer

Position Control: Optimization of height and angle, flame and hydride switch within one min

Background Correction

Background correction is available for both flame and graphite furnace methods. Deuterium lamp for flame. deuterium lamp or Zeeman for graphite furnace.

Correction Mode: Deuterium lamp, Zeeman Correction Capability:

a) Deuterium lamp: When background absorption approaches 1.0 Abs, the instrument is capable of a background correction of 60 times or more.

b) Zeeman : When background absorption approaches 2.0 Abs, the instrument is capable of a background correction of 100 times or more.

Data Processing:

Measurement Methods: Flame absorption, flame emission, graphite furnace, and hydride method.

Analysis Method: Linear fitting, nonlinear fitting, standard addition method.

Printing Output: Calibration curve, spectrum, analysis conditions, analysis parameters, and analysis results can be automatically stored and printed.

Dimensions: 1000 mm (L) x 610 mm (W) x 510 mm (H), 150 kg

Power: 220V / 50Hz, instrument 200W, graphite furnace 4 KW

HPLC Servicing, Validation, Trainings and Preventive Maintenance :

HPLC Servicing :HPLC Servicing : We have team of service engineers who can attend to any make of HPLC promptly @the most affordable cost.

Trainings :We also take up preventive Maintenance to reduce downtime of HPLC's Trainings.

AMC's/CMC :AMC's/CMC :We offer user training both in-House and at customer sites on HPLC principles, operations, trouble-shooting.

Validations :Validations :We have protocols for carrying out periodic Validations as per GLP/GMP/USFDA norms.

Instruments :Instruments :We offer instruments/Renting Services Modules like pumps,detector etc. on Rent.



About Analytical Technologies

Analytical Technologies is synonymous for offering technologies for doing analysis and is the Fastest Growing Global Brand having presence in at least 96 countries across the global. Analytical Technologies Limited is an ISO:9001 Certified Company engaged in Designing, Manufacturing, Marketing & providing Services for the Analytical, Chromatography, Spectroscopy, Bio Technology, Bio Medical, Clinical Diagnostics, Material Science & General Laboratory Instrumentation. Analytical Technologies, India has across the Country operations with at least 4 Regional Offices, 6 Branch Offices & Service Centers. Distributors & Channel partners worldwide.

Our Products & Technologies




UV/VIS
Spectro 2080+
Double Beam



Infra FTIR



Optima Gas
Chromatograph
3007

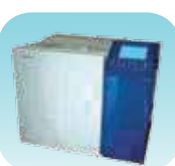

Optima Gas
Chromatograph
2979 Plus



Flash
Chromatograph



Atomic Absorption
Spectrophotometer



Liquid Particle
Counter



Optical Emission
Spectrophotometer



DSC/TGA



Semi Auto Bio
Chemistry Analyzer



HEMA 2062
Hematology
Analyzer



Micro Plate
Reader/Washer



URINOVA 2800
Urine Analyzer



Total Organic
Carbon 3800



Fully Automated
CLIA



NOVA-2100
Chemistry Analyzer



PCR/Gradient PCR/
RTPCR



TOC
Analyzer



Laser Particle
Size Analyzer



Ion Chromatograph



Water purification
system

Regulatory compliances



Corporate Social Responsibility

Analytical Foundation is a nonprofit organization (NGO) found for the purpose of:



1. Research & Innovation Scientist's awards/QC Professional Award : Quality life is possible by innovation only and the innovation is possible by research only, hence ANALYTICAL FOUNDATION is committed to identify such personalities for their contributions across various field of Science and Technology and awarding them yearly. To participate for award, send us your details of research / testing / publication at Info@analyticalfoundation.org

2. Improving quality of life by offering YOGA Training courses, Work shops/Seminars etc.

3. ANALYTICAL FOUNDATION aims to DETOXYFY human minds,souls and body by means of yoga, Meditation, Ayurveda, Health Care, Awards, Media, Events, Camps etc.

Reach us @



Note : Company reserves rights to add/delete/modify the contents / technical specifications of the catalogue without prior notice.

HPLC Solutions MultipleLabs Analytical Bio-Med Analytical Distributors Analytical Foundation (Trust)

Corporate & Regd. Office:
Analytical House, # E67 & E68,
Ravi Park, Vasna Road, Baroda,
Gujarat 390 015. INDIA

T: +91 265 2253620
+91 265 2252839
+91 265 2252370
F: +91 265 2254395

E: info@hplctechnologies.com
info@multiplelabs.com
info@analyticalgroup.net
info@analyticalbiomed.com

W. www.analyticalgroup.net
www.hplctechnologies.com
www.multiplelabs.com
www.ais-india.com

Sales & Support Offices:
across the country :
Distributors & Channel
partners World Wide