

CSP - 3800SP

Cross Section Polisher



EPCC / PRODUCTS / APPLICATION / SOFTWARE / ACCESSORIES / CONSUMABLES / SERVICES

Analytical Technologies Limited

An ISO 9001 Certified Company


www.analyticalgroup.net

- The CSP-3800SP is an advanced sample preparation tool that etches a cross section of a sample using an argon ion beam. This process avoids physical deformation and structural damage, without requiring complicated chemical processes. In addition, the system simplifies cross-sectional analysis of the sample by processing large areas from tens of μm to several mm.

►► FEATURES:

- A high etching rate of 700 μm per hour (based on Si, 8 kV)
- Ability to save/load recipes that are frequently used
- Step by step recipes with automatic execution function
- Easy to load sample using smart sample holder
- Real-time observation of ion beam status and etching status through chamber camera
- Convenient operation - intuitive GUI and easy touch screen
- Minimizes thermal damage with the ion beam Auto On/Off function
- Fast and convenient sample alignment with ion beam using built-in digital microscope
- Provided with noise, vibration, oil-free diaphragm pump
- Flat milling function provided for plane etching of a large areas

►► PRINCIPLES OF CROSS SECTION POLISHER:

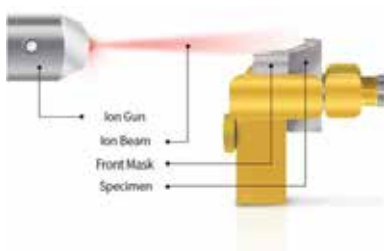


When a voltage is applied to the ion gun and argon gas is injected, plasma is generated and an ion beam is directed at the sample by an acceleration voltage to begin the etching process. If the sample is located behind a metal mask and the ion beam is directed at the metal mask and sample, the shielding effect of the metal mask minimizes beam damage to obtain clean cross-sectional etching results.

►► MECHANICAL POLISHING VS CP POLISHING:

If you polish using a mechanical polishing device, it is difficult to check the exact state of the cross section due to physical damage and contamination, but when performing cross-section processing with CP using an ion beam, you can observe the micro-surface structure of the sample without structural damage and contamination.




Mechanical Polishing Only

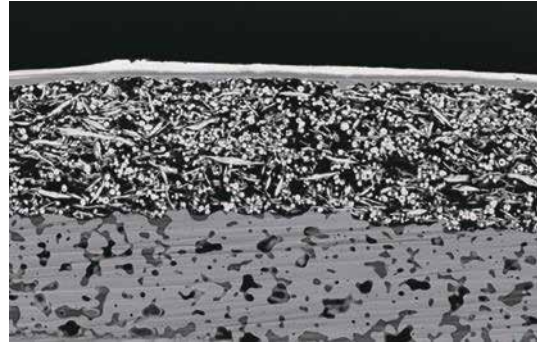


CP Polishing

►► FIB POLISHING VS CP POLISHING:

When etching a single side of the same sample with the CSP-3800SP, you can achieve excellent time and cost savings by etching a much wider cross section in a shorter time, compared to focused ion beam (FIB). Following images are the results of the same specimen milled in two different ways, FIB and CP for the same period of time.

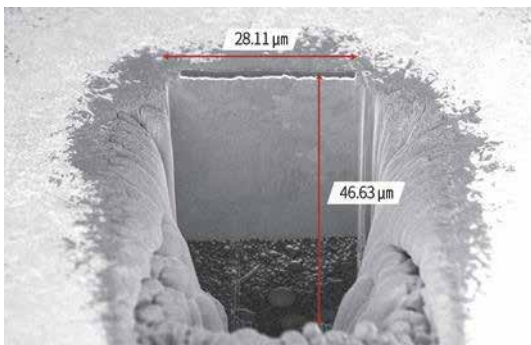


FIB Polishing

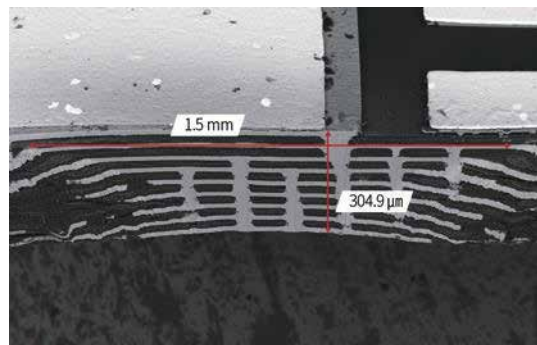

CP Polishing

►► CP PROCESS:

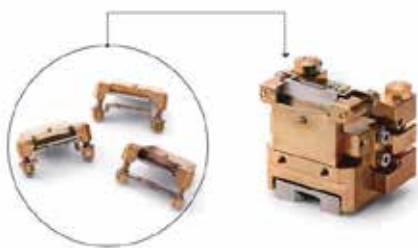


1. Select a holder that fits the size of the sample and mount it on the sample stand.



2. Adjust the height of the sample by selecting the sample height jig.
(Provide with 30, 50, 80 μm jig)




3. Fix the sample stand to the digital microscope.

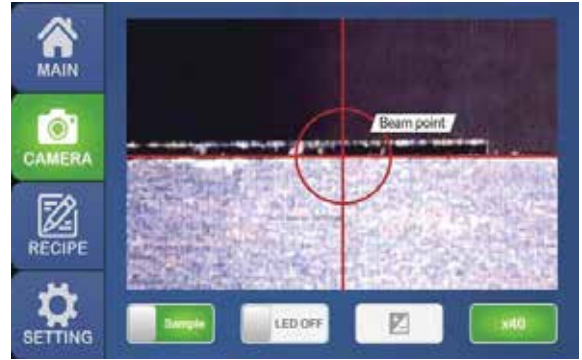

4. Select Sample Camera on the screen and place the desired etching position in the middle of the cross line.



5. Open the chamber door and mount the sample stand on the stage.

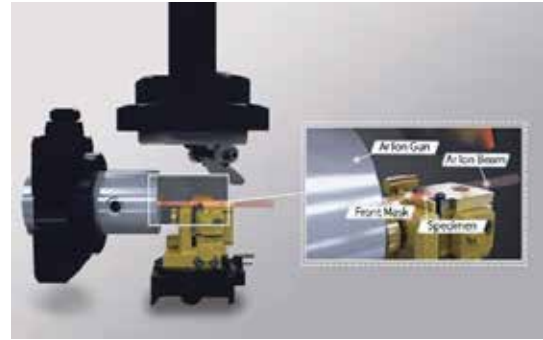

6. After entering the desired settings on the screen, press Run to begin etching.

►► INTUITIVE GUI:



1 Camera
: Digital Microscope / Chamber Camera


2 Auto Beam On/Off Mode

3 Step-by-step Mode

4 Flat Milling Mode

▶▶ DIGITAL MICROSCOPE:

Using a digital microscope, the position of the ion beam and the position of the sample can be easily aligned through the screen.

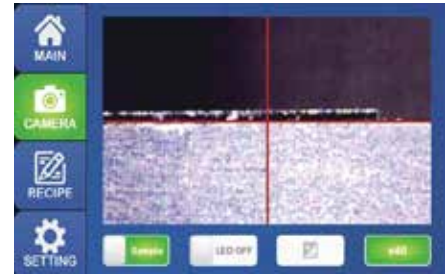

▶▶ CHAMBER CAMERA:


You can check the status of the ion beam through the camera inside the chamber and check the degree of cross-sectional etching in real time.



▶▶ AUTO BEAM ON / OFF MODE:

A feature designed to minimize heat damage caused by the ion beam by turning the beam on and off according to the set Ion Beam On/Off timer. It is useful for obtaining accurate cross-sectional conditions when etching heat-sensitive samples such as polymers and paper.

Paper >>

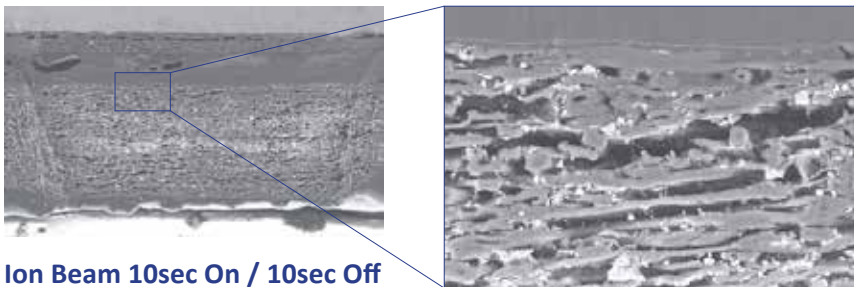


▶▶ AUTO BEAM ON / OFF MODE:

Frequently used etching conditions can be stored in the recipe list to easily apply setting values whenever necessary. In addition, a step-by-step mode is also available to store several recipes and automatically execute them to etch samples.





Before using STEP BY STEP MODE
(5kV 1hour)

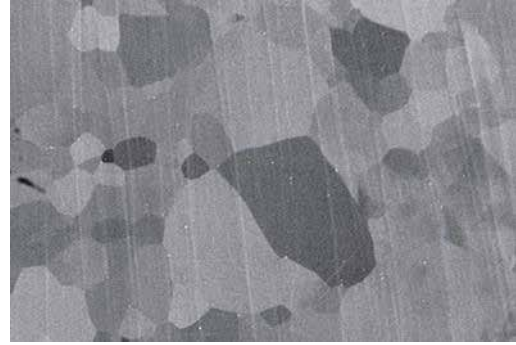


After using STEP BY STEP MODE
(5kV 1hour + 3kV 30min)

►► FLAT MILLING:

The CSP-3800SP can process samples in a plane using a dedicated holder. When a sample is mounted in the dedicated holder and the flat milling function is used, several mm² areas are etched by the ion beam based on the rotation center axis. At this point, since the polishing speed, area, and depth vary depending on the angle of incidence at which the ion beam hits the sample surface, uniform surface polishing should be induced by rotating and adjusting the angle of the sample. As the ion beam irradiates a larger area to etch an oxide layer or foreign substance, it is useful for pre-treatment of a large area of the sample.

MgO >>

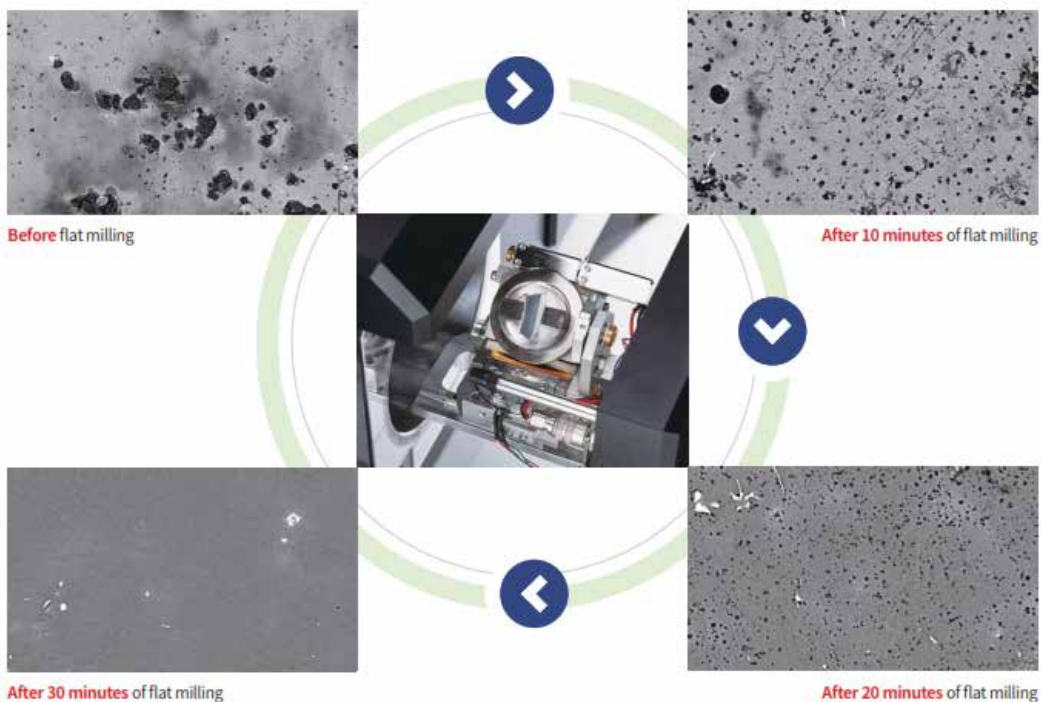


	Before milling	After milling
Element	Weight (%)	
C	12.31	6.27
O	29.44	0.59
Mg	56.42	92.25
Al	1.20	-
Mn	0.10	0.21
Zn	0.54	0.68
Total	100.00	100.00



►► CRYSTALLOGRAPHY OF METALLIC MATERIAL :

Precise surface pre-treatment is a very important factor for accurate grain and grain boundary analysis of metal samples. Ion beam milling is an ideal method to prepare samples prior to studying the mechanical and electrical properties of the sample through EBSD analysis.


Cu >>

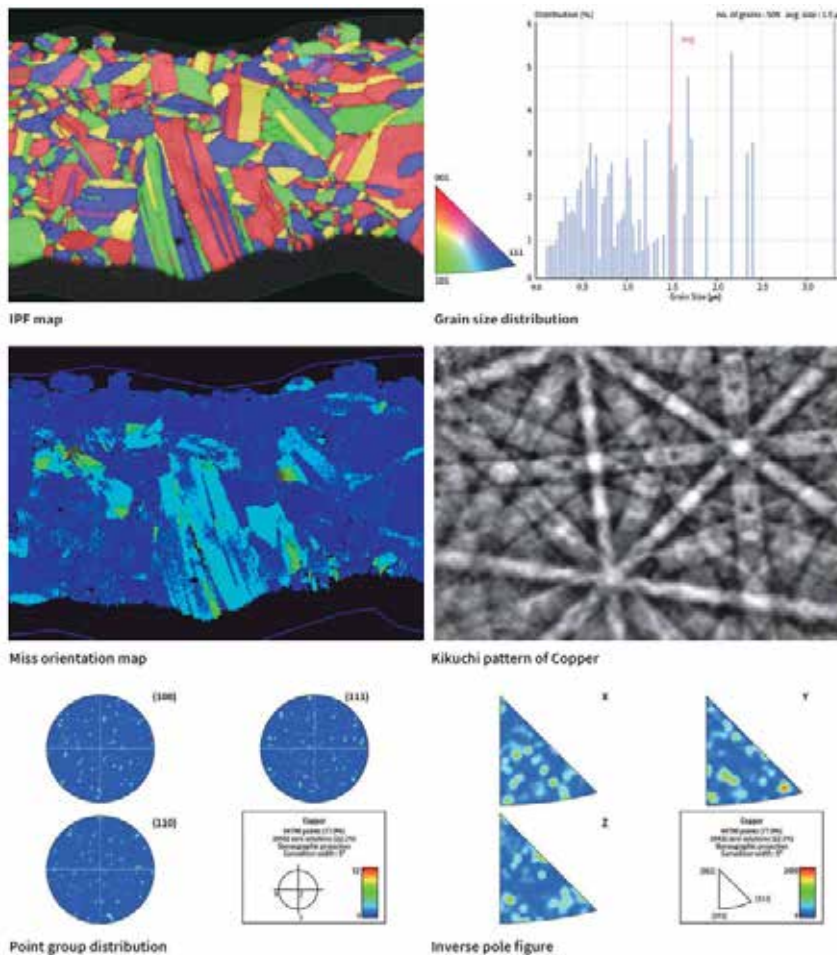



►► SEMI-CONDUCTOR

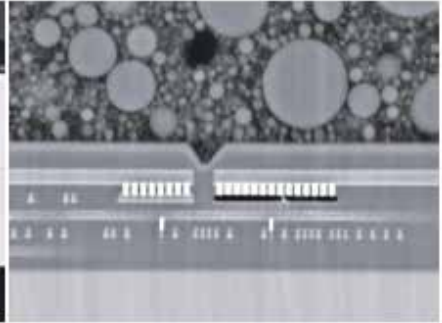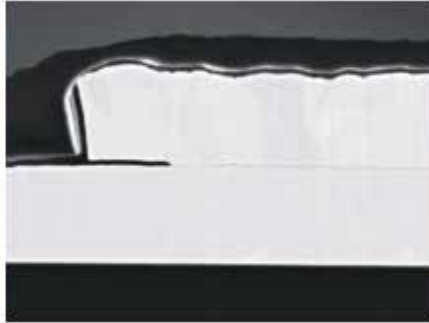

►► POWDER

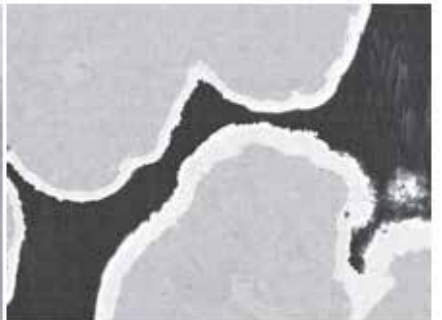



►► SOLAR CELL

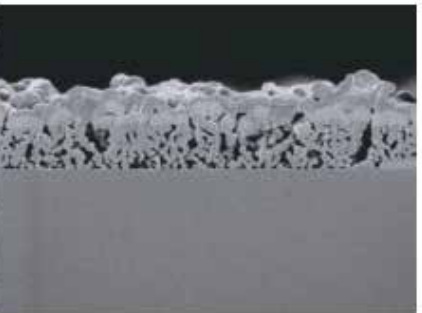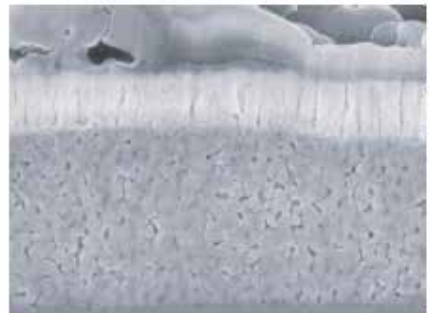


►► BATTERY




►► FIBER

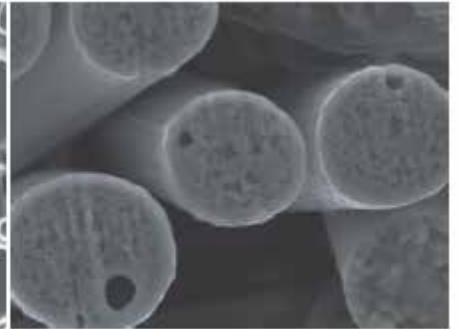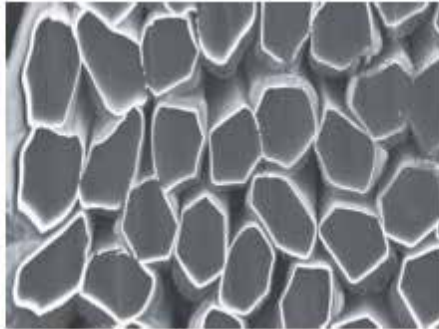


►► MINERALS

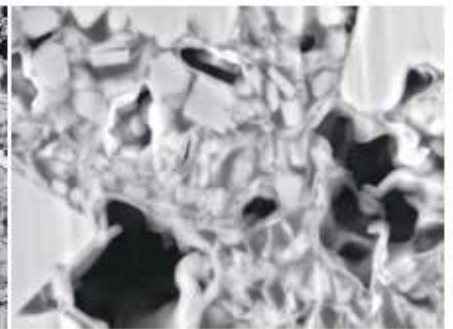

►► ALLOYS

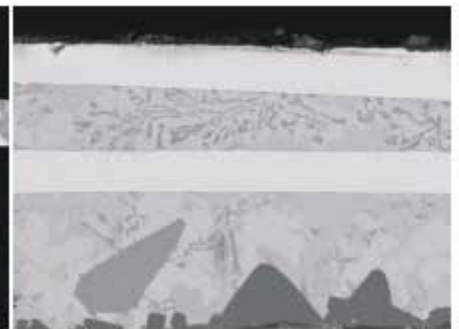


►► FILMS

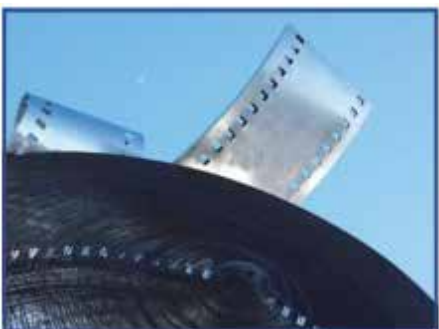

►► CRYSTALLOGRAPHY OF METALLIC MATERIAL :

After removing the oxide layer and foreign substances in the sample cross section through ion milling, EPMA component analysis for more accurate qualitative and quantitative analysis is possible. The following is a mineral sample, which was etched with CP to remove foreign substances and oxide layers on the surface before analyzing the elemental distribution through EPMA analysis.


Minerals >>

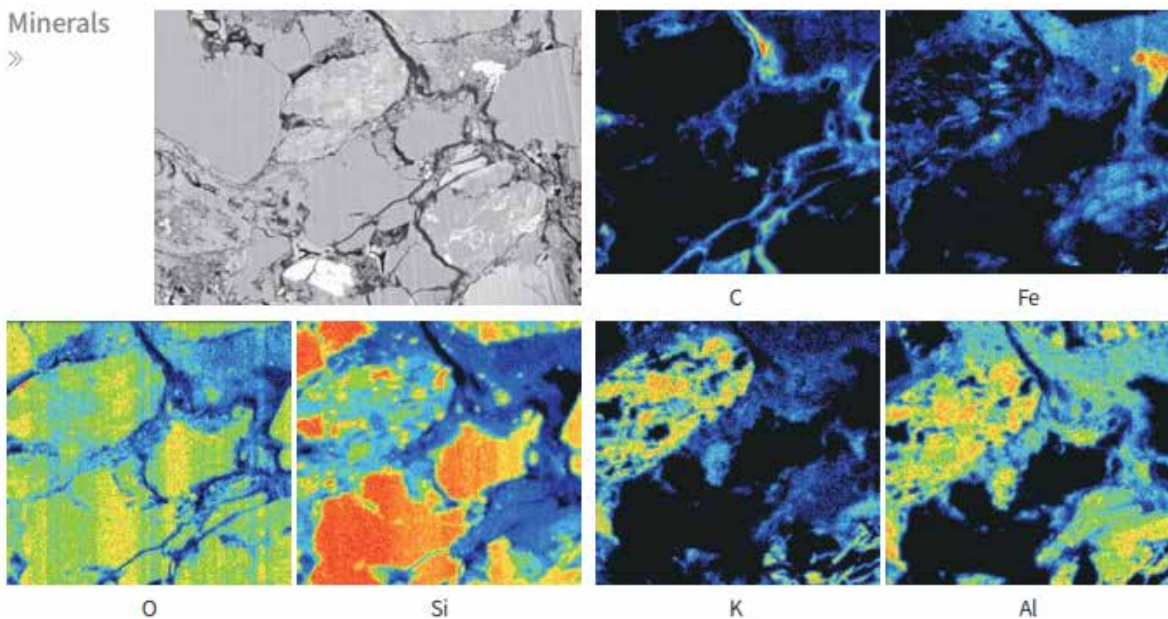

►► EDS :

By etching a sample with an ion beam to minimize damage to the cross-sectional structure and analyzing the components with EDS, the exact distribution of surface components can be observed without structural distortion of the sample.

MLCC
 (Multi layer ceramic capacitor)

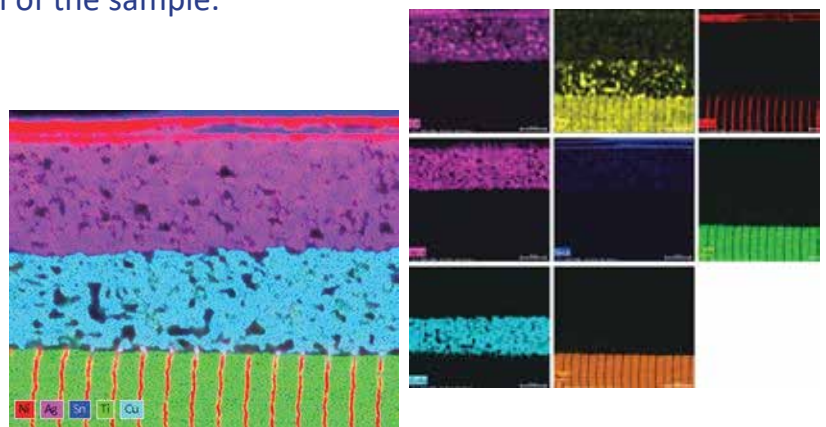

►► SPECIFICATIONS:

Accelerating voltage :	2 to 8kV
Milling rate :	700 μ m/h (at 8kV on Si wafer)
Sample stage swing angle :	$\pm 35^{\circ}$
Maximum sample size :	20(W) \times 10(L) \times 5.5(T)mm 16(W) \times 10(D) \times 9.5(H)mm
Specimen movement range :	X axis movement : ± 1.5 mm / Y axis movement : ± 2 mm
Flat milling stage tilt angle range :	40° to 80°
Sample size for flat milling :	$\varnothing 30 \times 11.4$ (H)mm
Operation :	7 inch touch panel
Digital Microscope for sample positioning :	Mag. x5, x10, x20, x40
Chamber camera for monitoring :	Mag. x5, x10, x20, x40 Brightness adjustable in 4 steps Ion beam observation mode (LED Off)
Gas for Ion :	Argon gas (99.999%)
Gas pressure :	0.1 Mpa (14.5psi)
Gas flow control :	Mass Flow Control
Vacuum systems :	Turbo pump, Diaphragm pump
Dimension :	607(W) \times 472(D) \times 277.5(430.5)(H) mm
Weight :	Main system 36kg / Diaphragm pump 6.5kg
Features :	Auto Beam On/Off mode Step by step mode

HPLC Servicing, Validation, Trainings and Preventive Maintenance :

HPLC Servicing :HPLC Servicing : We have team of service engineers who can attend to any make of HPLC promptly @the most affordable cost.

Trainings :We also take up preventive Maintenance to reduce downtime of HPLC's Trainings.

AMC's/CMC :AMC's/CMC :We offer user training both in-House and at customer sites on HPLC principles, operations, trouble-shooting.

Validations :Validations :We have protocols for carrying out periodic Validations as per GLP/GMP/USFDA norms.

Instruments :Instruments :We offer instruments/Renting Services Modules like pumps,detector etc. on Rent.



About Analytical Technologies

Analytical Technologies is synonymous for offering technologies for doing analysis and is the Fastest Growing Global Brand having presence in at least 96 countries across the global. Analytical Technologies Limited is an ISO:9001 Certified Company engaged in Designing, Manufacturing, Marketing & providing Services for the Analytical, Chromatography, Spectroscopy, Bio Technology, Bio Medical, Clinical Diagnostics, Material Science & General Laboratory Instrumentation. Analytical Technologies, India has across the Country operations with at least 4 Regional Offices, 6 Branch Offices & Service Centers. Distributors & Channel partners worldwide.

Our Products & Technologies



UV/VIS
Spectro 2080+
Double Beam



Infra FTIR



Optima Gas
Chromatograph
3007



Optima Gas
Chromatograph
2979 Plus



Flash
Chromatograph



Atomic Absorption
Spectrophotometer



Liquid Partical
Counter



Optical Emission
Spectrophotometer



DSC/TGA



Semi Auto Bio
Chemistry Analyzer



HEMA 2062
Hematology
Analyzer



Micro Plate
Reader/Washer



URINOVA 2800
Urine Analyzer



Total Organic
Carbon 3800



Fully Automated
CLIA



NOVA-2100
Chemistry Analyzer



PCR/Gradient PCR/
RTPCR



TOC
Analyzer



Laser Particle
Size Analyzer



Ion Chromatograph



Water purification
system

Regulatory compliances



Corporate Social Responsibility

Analytical Foundation is a nonprofit organization (NGO) found for the purpose of:



Analytical
Foundation

1. Research & Innovation Scientist's awards/QC Professional Award : Quality life is possible by innovation only and the innovation is possible by research only, hence ANALYTICAL FOUNDATION is committed to identify such personalities for their contributions across various field of Science and Technology and awarding them yearly. To participate for award, send us your details of research / testing / publication at Info@analyticalfoundation.org

2. Improving quality of life by offering YOGA Training courses, Work shops/Seminars etc.

3. ANALYTICAL FOUNDATION aims to DETOXYFY human minds,souls and body by means of yoga, Meditation, Ayurveda, Health Care, Awards, Media, Events, Camps etc.

Reach us @



 **Analytical**[®]
Technologies Limited

HPLC Solutions MultipleLabs Analytical Bio-Med Analytical Distributors Analytical Foundation (Trust)

Corporate & Regd. Office:
Analytical House, # E67 & E68,
Ravi Park, Vasna Road, Baroda,
Gujarat 390 015. INDIA

T: +91 265 2253620
+91 265 2252839
+91 265 2252370
F: +91 265 2254395

E: info@hplctechnologies.com
info@multiplelabs.com
info@analyticalgroup.net
info@analyticalbiomed.com

W. www.analyticalgroup.net
www.hplctechnologies.com
www.multiplelabs.com
www.ais-india.com

Sales & Support Offices:
across the country :
Distributors & Channel
partners World Wide

Note : Company reserves rights to add/delete/modify the contents / technical specifications of the catalogue without prior notice.