

OM - 3216 Osmometer



EPCC / PRODUCTS / APPLICATION / SOFTWARE / ACCESSORIES / CONSUMABLES / SERVICES

Analytical Technologies Limited

An ISO 9001 Certified Company

www.analyticalgroup.net



►► Introduction

- Sample volume 10 to 25 µl · Measuring time approx. 1,3 minutes
- Simple operation with touch-display and Icon-style
- Air cooling Clear sample identification with built in scanner and time clock Single-use plastic sample tubes
- 3-point variable calibration Built in printer · Multi level user access

►► Measuring Principle

The freezing point of aqueous solutions is measured. The depression of freezing point compared to pure water is direct related to the osmotic concentration. Pure water freezes at 0°C, an aqueous solution with an osmolality of 1 osmol / kg water at -1.858°C. 1 Mol of a substance dissolved in 1 kg of water gives a solution with an osmotic concentration of 1 osmol / kg water only if it is an ideal solution and if the substance does not dissociate.

Typical cooling curves of water and aqueous solutions

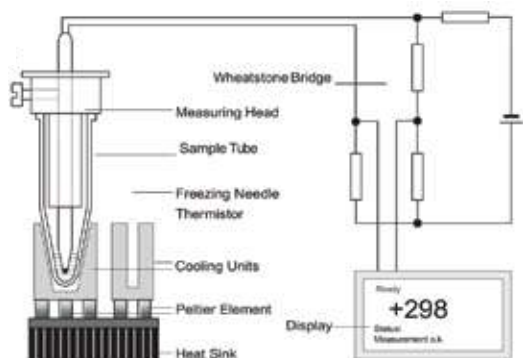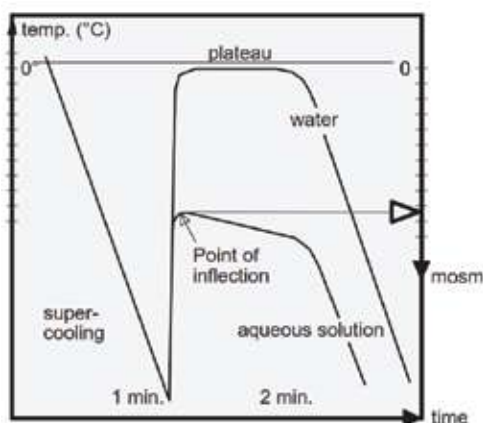

Schematic diagram

►► **Function and Description**

The sample (serum, blood, urine or any other aqueous solution) is filled in a plastic sample tube and cooled via a Peltierelement. These are semiconductor devices, which become cold on one side and warm on the other when electrical current flows through. The cold side cools the sample whilst the heat from the warm side dissipates into the air by convection.

The Osmometer transports the warmth via cooling surfaces into the atmosphere, an extra water supply is not necessary. Operation of the instrument requires only an electrical socket. The temperature on the cold side is kept electronically constant.

During the measuring process the temperature of the sample is measured by a thermistor (a temperature dependent resistor). This is part of the measuring head onto which the tube is placed.

The measuring head is attached to guide rods which protect it from accidental damage.

At a defined supercooling the freezing process is started by lowering a needle with ice crystals into the sample tube. The freezing point of the sample is reached.

The method of initiating the freezing process is important for the reproducibility of measurements. Dipping a needle with ice crystals into the sample gives more exact results than stirring with a wire which is constantly dipping into the sample. Because of linear correlation between osmolality and freezing point the measurement of freezing point is a determination of osmolality. The results are displayed as mosm / kg H₂O.

►► **Technical Data**

Sample volume: 10 to 50 µl

Measuring time: approx. 1.3 minutes (25 µl)

Reproducibility: ± 0,5% (50 µl), ± 1 % (10 to 25 µl)

Measurement range: 0 ... 2500 mosm / kg H₂O

Measurement display changeable (mosm / m°C/%0,9NaCl)

Measurement memory: 200 measurements with sample numbers

Data interfaces: USB/RS 232 for PC and RS232 for handscanner

Power supply: 100-230V AC approx. 65VA

Dimensions: Width 190 mm, height 278 mm, depth 216 mm

Weight: approx. 4.3 kg

The Instrument is CE-labelled

►► Accessories included

50 plastic tubes, 2x10 vials of standard solution 300 and 900 mosm / kg H₂O, Spare thermistor, spare needle, spare fuses, Cleaning fluid and little funnel for cleaning needle hole, USB/RS232 cables and data transmitting program, operating instructions.

►► Calibration

The zero point is calibrated with distilled water and a standard point with the NaCl-solution of 300 mosm / kg H₂O supplied. Occasional checks shows that the calibration is very stable. An additional 900 mosm-solution can be used to calibrate, whereby the linearity is raised by higher concentrations.

The second and third calibration point could be changed by user to meet other standard concentrations.

►► Special Features and Advantages

- Touch operation with Icon-style on a wide black & white graphic LC display.
- Selectable languages.
- Multi level user access with passwords and usernames.
- Single use plastic tubes.
- Blocking of measuring tube or needle by freezing in stand-by mode is counteracted by regular automatically defrosting.
- Needle can be rinsed with water or disinfection solution.
- Built in thermoprinter and scanner for sample numbers.
- Integrated real time clock to sort measurements and calibrations.
- Reduced operating costs as no extra water supply is necessary.
- Wide input voltage range 100-230V AC suitable for all countries.

►► Operating of the Instrument

- Switch on instrument, wait for ready display (approx.3 minutes).
- Place 25 µl sample in the sample tube ($\pm 10\%$).
- Place the sample tube onto the measuring head.
- Lower the measuring head. The sample tube is dipped into the cooling slot.
- Enter sample number using the built in scanner or touch-display. These entries can also be made after measurement.

- Superercooling is reached after 1 minute and the needle with the ice crystals is dipped into the sample.
- As freezing point is reached, the display shows the measurement value in mosm / kg and the instrument gives an audible tone. The measurement value is stored in the memory.
- Slide measuring head from cooling slot and remove the sample tube. Result will be printed with the built in printer.
- Wipe the thermistor with a soft tissue.

►► **Extra Accessories**

- Analogue output for XY-plotter
- Calibration solution 300 and 900 mosm / kg H₂O

HPLC Servicing, Validation, Trainings and Preventive Maintenance :

HPLC Servicing :HPLC Servicing : We have team of service engineers who can attend to any make of HPLC promptly @the most affordable cost.

Trainings :We also take up preventive Maintenance to reduce downtime of HPLC's Trainings.

AMC's/CMC :AMC's/CMC :We offer user training both in-House and at customer sites on HPLC principles, operations, trouble-shooting.

Validations :Validations :We have protocols for carrying out periodic Validations as per GLP/GMP/USFDA norms.

Instruments :Instruments :We offer instruments/Renting Services Modules like pumps,detector etc. on Rent.



About Analytical Technologies

Analytical Technologies is synonymous for offering technologies for doing analysis and is the Fastest Growing Global Brand having presence in at least 96 countries across the global. Analytical Technologies Limited is an ISO:9001 Certified Company engaged in Designing, Manufacturing, Marketing & providing Services for the Analytical, Chromatography, Spectroscopy, Bio Technology, Bio Medical, Clinical Diagnostics, Material Science & General Laboratory Instrumentation. Analytical Technologies, India has across the Country operations with at least 4 Regional Offices, 6 Branch Offices & Service Centers. Distributors & Channel partners worldwide.

Our Products & Technologies



UV/VIS
Spectro 2080+
Double Beam



Infra FTIR



Optima Gas
Chromatograph
3007



Optima Gas
Chromatograph
2979 Plus



Flash
Chromatograph



Atomic Absorption
Spectrophotometer



Liquid Particle
Counter



Optical Emission
Spectrophotometer



DSC/TGA



Semi Auto Bio
Chemistry Analyzer



HEMA 2062
Hematology
Analyzer



Micro Plate
Reader/Washer



URINOVA 2800
Urine Analyzer



Total Organic
Carbon 3800



Fully Automated
CLIA



NOVA-2100
Chemistry Analyzer



PCR/Gradient PCR/
RTPCR



TOC
Analyzer



Laser Particle
Size Analyzer



Ion Chromatograph



Water purification
system

Regulatory compliances



Corporate Social Responsibility

Analytical Foundation is a nonprofit organization (NGO) found for the purpose of:



Analytical
Foundation

1. Research & Innovation Scientist's awards/QC Professional Award : Quality life is possible by innovation only and the innovation is possible by research only, hence ANALYTICAL FOUNDATION is committed to identify such personalities for their contributions across various field of Science and Technology and awarding them yearly. To participate for award, send us your details of research / testing / publication at Info@analyticalfoundation.org

2. Improving quality of life by offering YOGA Training courses, Work shops/Seminars etc.

3. ANALYTICAL FOUNDATION aims to DETOXYFY human minds,souls and body by means of yoga, Meditation, Ayurveda, Health Care, Awards, Media, Events, Camps etc.



 **Analytical**®
Technologies Limited

HPLC Solutions MultipleLabs Analytical Bio-Med Analytical Distributors Analytical Foundation (Trust)

Corporate & Regd. Office:
Analytical House, # E67 & E68,
Ravi Park, Vasna Road, Baroda,
Gujarat 390 015. INDIA

T: +91 265 2253620
+91 265 2252839
+91 265 2252370
F: +91 265 2254395

E: info@hplctechnologies.com
info@multiplelabs.com
info@analyticalgroup.net

W. www.analyticalgroup.net
www.hplctechnologies.com
www.multiplelabs.com

Sales & Support Offices:
across the country :
Distributors & Channel
partners World Wide